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
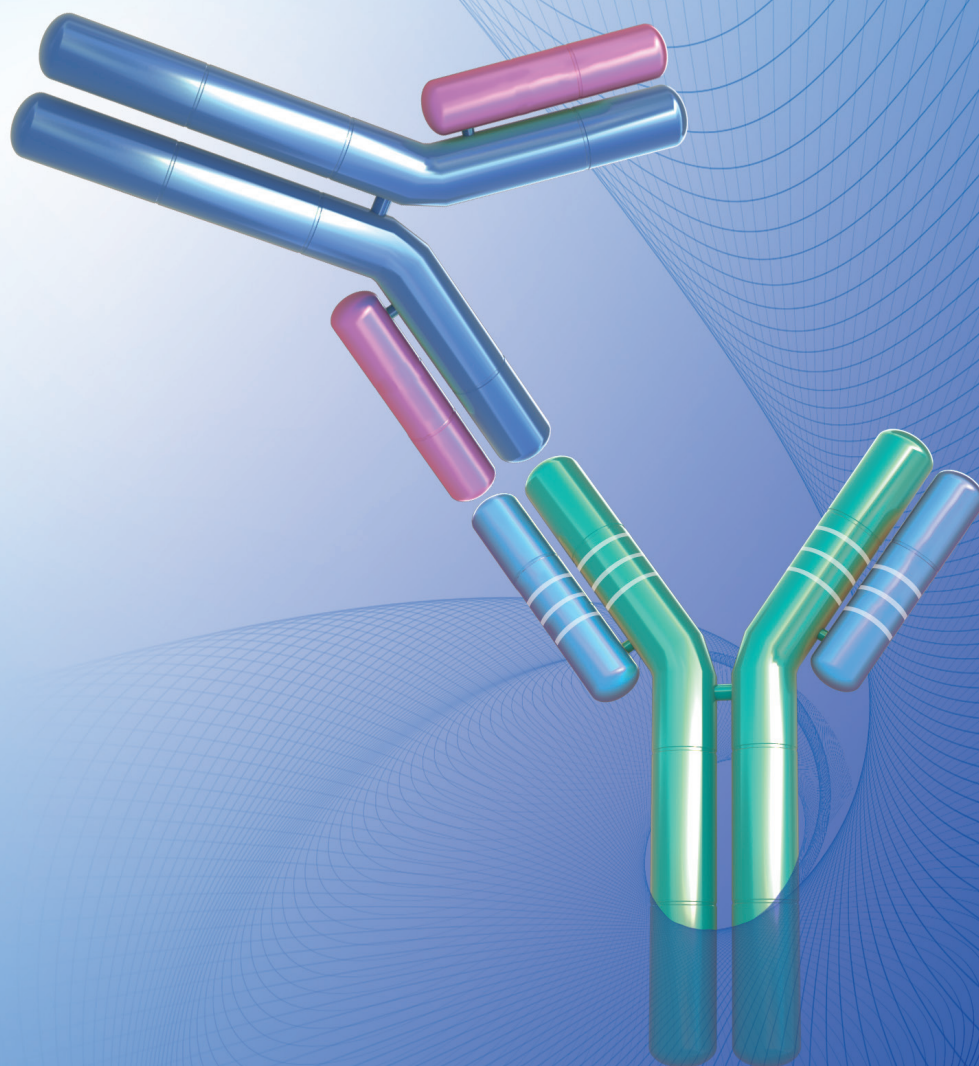
Figure 1. Example of Bridging ELISA-based PK assay

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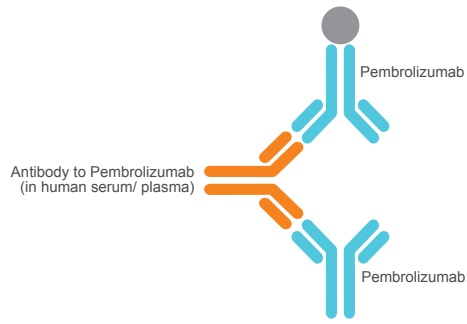Figure 2. Example of Bridging ELISA-based ADA assay

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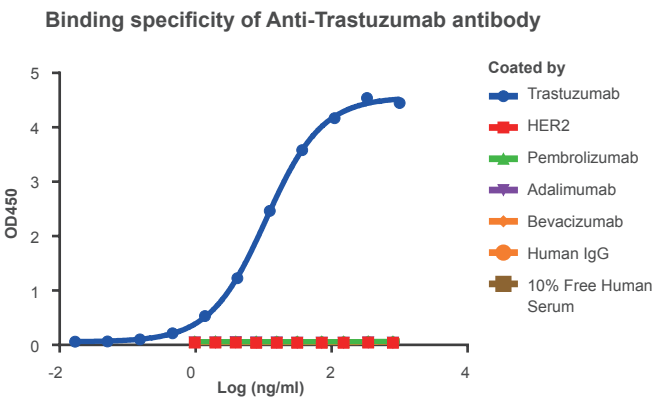
Figure 3. High specificity of Anti-Trastuzumab antibody to Trastuzumab.

The binding specificity of Anti-Trastuzumab Antibody (15H2), mAb, Mouse (A02033) to Trastuzumab was studied using indirect ELISA. No cross-reactivity was noted with human IgG, 10% free human serum, and other antibody drugs shown in the figure.
Note: 10% free human serum refers to T3/T4 Free human serum, lipid stripped.

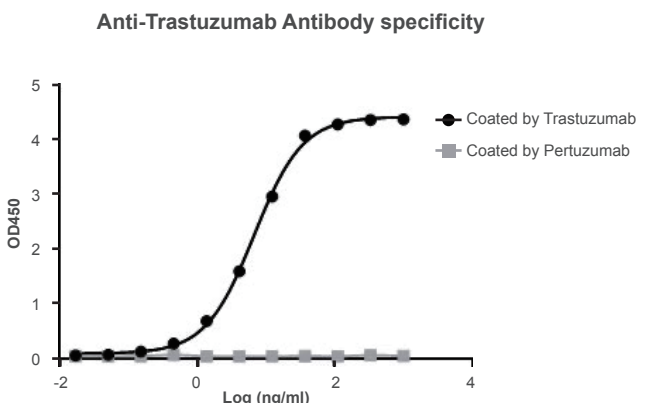
Figure 4. High specificity of Anti-Trastuzumab to Trastuzumab.

The binding specificity of Anti-Trastuzumab Antibody (15H2), mAb, Mouse (A02033) to Trastuzumab was studied using indirect ELISA. No cross-reactivity with Pertuzumab was observed.

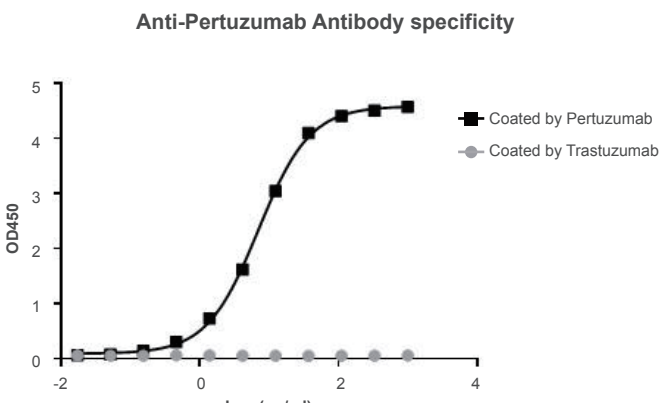
Figure 5. High specificity of Anti-Pertuzumab to Pertuzumab.

The binding specificity of Anti-Pertuzumab Antibody (1A3), mAb, Mouse (A02150) to Pertuzumab was studied using indirect ELISA. No cross-reactivity with Trastuzumab was observed.

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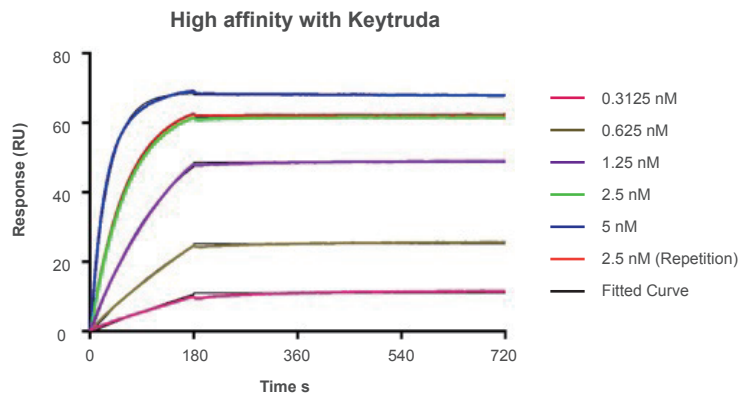
Figure 6. High affinity of Anti-Pembrolizumab antibody with Keytruda.

The affinity of MonoRab™ Anti-Pembrolizumab Antibody (90G12F8), mAb, Rabbit (Cat.No. A01853) towards Keytruda was measured by SPR method; $K_d < 10^{-12}$ M.

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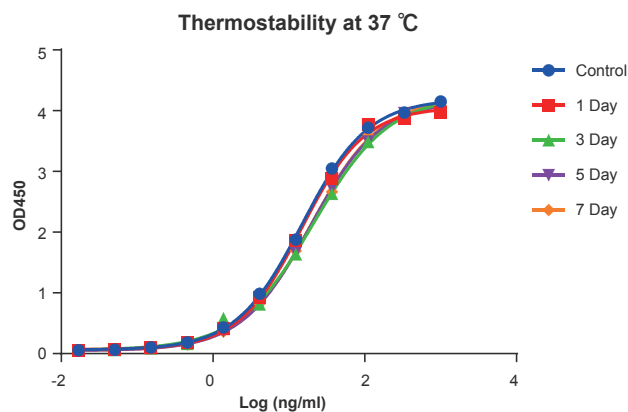
Figure 7. High thermal stability of A02033.

Anti-Trastuzumab Antibody (15H2), mAb, Mouse (A02033, 0.5 mg/ml) was aliquoted and placed at 37°C for different time periods (1, 3, 5, 7 days), along with the control placed at 4°C. No significant loss of binding activity was observed. Trastuzumab (1 µg/ml) was used as coating antigen.

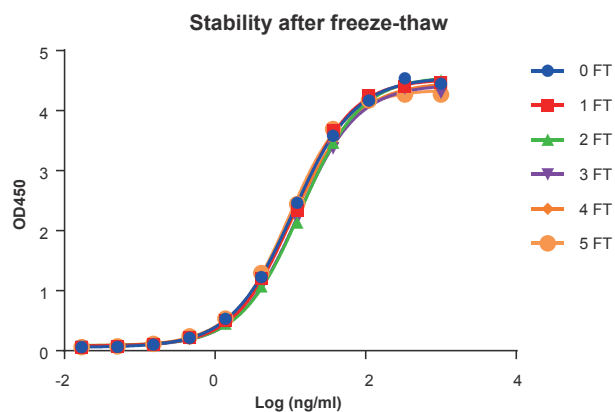
Figure 8. High freeze-thaw stability of A02033.

Anti-Trastuzumab Antibody (15H2), mAb, Mouse (A02033, 0.5 mg/ml) were subjected to the indicated number of freeze-thaw cycles (FT). No significant loss of activity was observed.

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